METTLER TOLEDO

O₂ ISM Qualification Kit

**Instruction manual
Bedienungsanleitung
Instructions d'utilisation**

O₂ ISM Qualification Kit
52 300 440

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Instruction manual

Contents

	Page
1 Safety.....	5
2 Product Description	7
3 Operation	9
3.1 Connecting the transmitter to the O ₂ ISM Qualification Kit	9
4 Product Specification	12
4.1 Simulator, Pre-registered Values	12
4.2 Sensor Wear Data.....	13
4.3 Warnings and Alarms	15
5 Installation	16
5.1 Dimensions	16
6 Ordering Information	17
7 Certificates	18
8 Protocol.....	19

1 Safety

This instruction manual lists the main points of how to use the **O₂ ISM Qualification Kit**, from METTLER TOLEDO for best results and in the intended manner.

This **instruction manual**, and the **safety requirements** in particular, **must be observed** by the **operator and all other persons** working with this equipment.

The instruction manual must always be stored close at hand, in a place accessible to all people working with an ISM O₂ sensor and the O₂ ISM Qualification Kit.

Warning notices and symbols

This instruction manual identifies safety instructions and additional information by means of the following symbols:

This symbol draws attention to **safety instructions and warnings of potential danger** which, if neglected, could result in injury to persons and/or damage to property.

This symbol identifies **additional information and instructions** which, if neglected, could lead to defects, inefficient operation and possible loss of product.

Intended use

The O₂ ISM Qualification Kit is intended **solely for validating and testing of the ISM transmitter family**, respectively, as described in this instruction manual.

The O₂ ISM Qualification Kit is intended **for simulating the functions of METTLER TOLEDO ISM sensors**, as described in this instruction manual.

Any other use, or use not mentioned here, that is incompatible with the technical specifications is deemed inappropriate. The operator is solely responsible for any damage arising from such use.

Other prerequisites for appropriate use include:

- observing the instructions, notes and requirements set out in this instruction manual
- correct care and maintenance of the unit, according to the instruction manual
- observing the prescribed environmental and operational conditions
- compliance with local statutes

2 Product Description

The O₂ ISM Qualification Kit is a set of 5 different sensor plug-in heads and allows the simulation of the following sensor types with predefined values and errors (not changeable by the user) to assure safe use as qualification tool:

- InPro 6850 i
- InPro 6900 i
- InPro 6950 i

Each sensor plug-in head corresponds to one of the mentioned METTLER TOLEDO sensors and delivers the complete data of an ISM sensor.

The O₂ ISM Qualification Kit has basically 2 functions:

- Tool for the qualification of measuring systems.
- Service tool for quick checks and loop trouble-shooting.

The data indicated (primary/secondary) and transmitted are as follow:

- Main measured value
- Temperature
- Error message according to indication.

All standard process and wear data of the sensor are also available in each plug-in head and are also sent to the transmitter or the PLC as well as to the iSense software.

3 Operation

3.1 Connecting the transmitter to the O₂ ISM Qualification Kit

To qualify your measuring systems proceed as follows:

Please refer to the transmitter instruction manual for the exact description of the transmitter operation.

Step 1: Disconnect the ISM sensor from the measuring point.

Step 2: Connect one of the O₂ ISM Qualification Kit plug-in heads to the transmitter via the standard connecting cable.

Step 3: The preregistered values then are displayed on the transmitter.

Step 4: Check if the displayed values by the transmitter or the PLC correspond to the predefined fixed values of the O₂ ISM Qualification Kit plug-in head.

See certification for exact values and tolerances.

Step 5: In case you are using the Error simulators, so check that the appropriate errors are correctly displayed by the transmitter or the PLC.

Step 6: Report results into the attached spreadsheet "Protocol for O₂ ISM Qualification Kit".

Please note: the O₂ ISM Qualification Kit simulates the behavior of a sensor at several temperatures. All commonly used tables programmed in each transmitter for solubility, salinity correction, etc. apply standardized values, e.g. temperature 25 °C (77 °F), pressure 1.013 bar (14.69 psi) and a relative humidity 100 %. The transmitter calculates back the value that the sensor would deliver under these reference conditions. This compensated value may differ from the direct current delivered by the simulator if for example the calibration pressure (manually tipped or automatically measured by some transmitter) differs from 1.013 bar (14.69 psi) or if the relative humidity typed in is different from 100 %.

Problem solving: If any of the tests described above failed, you should change the connecting cable and repeat the test. If the test still shows the same result, the transmitter or the cable is defective. If you are using a transmitter with current outputs, so ensure the outputs are correctly wired.

You must **never do a calibration** (the simulators are read-only).

You must **never increment autoclave cycles** (always set "NO" to cycle counter), this is only needed if autoclave has been set to a value higher than 0.

Different transmitters may display different values and errors.


Out of range behavior is also different depending on the transmitter in use thus **different values/readings may occur.**

Set the transmitter input filter to off (only for SIM O₂ TGL, SIM O₂ ppb TGL and SIM O₂ Trace TGL).

Do not use the O₂ ISM plug-in heads after the expiry date shown on the label of each plug-in head. Please send them back to METTLER TOLEDO for recertification.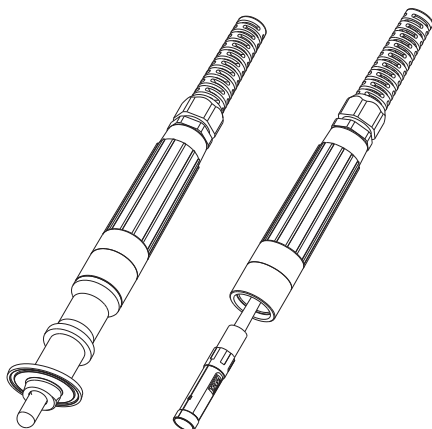

Measuring point with
METTLER TOLEDO Electrode

Connecting O₂ ISM
Qualifying plug-in head

Environment operating and storing conditions:

Permissible ambient temperature (operation):
–25 °C to 60 °C (–13 °F to 140 °F)

Permissible ambient temperature (storage/transport):
–25 °C to 60 °C (–13 °F to 140 °F)

4 Product Specification

4.1 Simulator, Pre-registered Values

Order No.	Product Designation	Legend	OXY Value	Temp.
Standard (InPro 6850i) SIM O₂ Kit			Order No. 52 300 416	
52 300 417	SIM O ₂ ZERO	SIM O ₂ ZERO	0%	25 °C
52 300 418	SIM O ₂ AIR	SIM O ₂ AIR	20.95 Vol% 100% (air) sat	40 °C
52 300 419	SIM O ₂ TGL	Toggle Mode 30 sec	0% & 100%	25 °C & 40 °C
52 300 420	SIM O ₂ ERR1	9999 % ... %sat	9999 % ... %sat	25 °C
52 300 421	SIM O ₂ ERR2	9999 % ... %sat	83 % ... %sat	90 °C
Standard (InPro 6900i) SIM O₂ ppb Kit			Order No. 52 300 422	
52 300 423	SIM O ₂ ppb ZERO	SIM O ₂ ppb ZERO	0%	5 °C
52 300 424	SIM O ₂ ppb AIR	SIM O ₂ ppb AIR	20.95 Vol% 100% (air) sat	25 °C
52 300 425	SIM O ₂ ppb TGL	Toggle Mode 30 sec	0% & 100%	5 °C & 25 °C
52 300 426	SIM O ₂ ppb ERR1	9999 % ... %sat	9999 % ... %sat	25 °C
52 300 427	SIM O ₂ ppb ERR2	9999 % ... %sat	83 % ... %sat	90 °C
Standard (InPro 6950i) SIM O₂ Trace Kit			Order No. 52 300 428	
52 300 429	SIM O ₂ Trace ZERO	SIM O ₂ Trace ZERO	0%	5 °C
52 300 430	SIM O ₂ Trace AIR	SIM O ₂ Trace AIR	20.95 Vol% 100% (air) sat	25 °C
52 300 431	SIM O ₂ Trace TGL	Toggle Mode 30 sec	0% & 100%	5 °C & 25 °C
52 300 432	SIM O ₂ Trace ERR1	9999 % ... %sat	9999 % ... %sat	25 °C
52 300 433	SIM O ₂ Trace ERR2	9999 % ... %sat	83 % ... %sat	90 °C

4.2 Sensor Wear Data

All O₂ SIM's except 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2), 52 300 433 (SIM O₂ TRACE ERR2)

Sensor Wear Monitor O₂	Transmitters M400/M420	% BAR
Time to maintenance	100 d	75 %
Life time indicator	2400 d	11 %
Adaptive Cal Timer	0 d	100 %
Sensor operating time	300 d	
Autoclaving cycles	1	
CIP cycles	2	
SIP cycles	3	
Max temperature	130/121 °C dd.mm.yyyy (date of production)	

For SIM 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2), 52 300 433 (SIM O₂ TRACE ERR2)

Sensor Wear Monitor O₂	Transmitters M400/M420	% BAR
Time to maintenance	0 d	100 %
Life time indicator	2300 d	15 %
Adaptive Cal Timer	0 d	100 %
Sensor operating time	300 d	
Autoclaving cycles	1	
CIP cycles	2	
SIP cycles	3	
Max temperature	130/121 °C dd.mm.yyyy (date of production)	

All O₂ SIM except 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2),
52 300 433 (SIM O₂ TRACE ERR2)

Sensor Wear Monitor O₂	Transmitter M700 %BAR
Membrane body wear	~100%
Inner body wear	~10%
Operating time	300 d
Network diagram sensor wear	50%

For SIM 52 300 421 (SIM O₂ ERR2), 52 300 427 (SIM O₂ PPB ERR2),
52 300 433 (SIM O₂ TRACE ERR2)

Sensor Wear Monitor O₂	Transmitter M700 %BAR
Membrane body wear	~100%
Inner body wear	~15%
Operating time	300 d
Network diagram sensor wear	25%

4.3 Warnings and Alarms


Warnings and Alarms	Transmitters M300/M400/M420
52 300 420 (SIM O ₂ ERR1)	Error O ₂ slope > -30 nA
52 300 426 (SIM O ₂ ppb ERR1)	Error O ₂ slope > -220 nA
52 300 432 (SIM O ₂ Trace ERR1)	Error O ₂ slope > -2000 nA
52 300 421 (SIM O ₂ ERR2)	Calibration required Maintenance required
52 300 427 (SIM O ₂ ppb ERR2)	Calibration required Maintenance required
52 300 433 (SIM O ₂ Trace ERR2)	Calibration required Maintenance required
Transmitter M700	
52 300 420 (SIM O ₂ ERR1)	Fail Measurement Range Warning Slope
52 300 426 (SIM O ₂ ppb ERR1)	Fail Measurement Range Warning Slope
52 300 432 (SIM O ₂ Trace ERR1)	Fail Measurement Range Warning Slope
52 300 421 (SIM O ₂ ERR2)	Fail Temperature Range Warning Sensor wear Warning Adjustment required
52 300 427 (SIM O ₂ ppb ERR2)	Fail Temperature Range Warning Sensor wear Warning Adjustment required
52 300 433 (SIM O ₂ Trace ERR2)	Fail Temperature Range Warning Sensor wear Warning Adjustment required

Please refer to your transmitter instruction manual for specific instructions.


5 Installation

5.1 Dimensions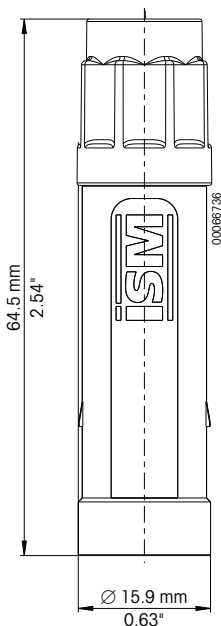

Markings

The markings labeled on each O₂ ISM plug-in head state the following information:

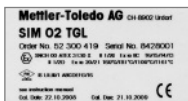6 Ordering Information

Order No.	Designation
52 300 410	SIM pH Kit
52 300 411	SIM pH4
52 300 412	SIM pH7
52 300 413	SIM pH TGL
52 300 414	SIM pH ERR1
52 300 415	SIM pH ERR2
52 300 416	SIM O₂ Kit
52 300 417	SIM O ₂ ZERO
52 300 418	SIM O ₂ AIR
52 300 419	SIM O ₂ TGL
52 300 420	SIM O ₂ ERR1
52 300 421	SIM O ₂ ERR2
52 300 422	SIM O₂ ppb Kit
52 300 423	SIM O ₂ ppb ZERO
52 300 424	SIM O ₂ ppb AIR
52 300 425	SIM O ₂ ppb TGL
52 300 426	SIM O ₂ ppb ERR1
52 300 427	SIM O ₂ ppb ERR2
52 300 428	SIM O₂ Trace Kit
52 300 429	SIM O ₂ Trace ZERO
52 300 430	SIM O ₂ Trace AIR
52 300 431	SIM O ₂ Trace TGL
52 300 432	SIM O ₂ Trace ERR1
52 300 433	SIM O ₂ Trace ERR2

The single simulators are not available for sale.

7 Certificates

ATEX/FM certification pending.

8 Protocol

Protocol for O₂ ISM Qualification Kit

Facility Name _____

Instrument Manufacturer _____

Date of Purchase _____

Transmitter Model _____

Transmitter Serial Nr _____

Simulator P/N	Simulator S/N	Designation	Valid until	Parameter/s (check all that apply for this instrument)

Date/Time	Acceptable? ⇒ Yes	Acceptable? ⇒ No	Name/Signature

Protocol for O₂ ISM Qualification Kit © Mettler-Toledo AG, MarCom, Im Hackacker 15, CH-8902 Urdorf. 03/2009 1

A copy of this form in A4 format is delivered with the O₂ ISM Qualification Kit.

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Protocole pour set de qualification ISM O₂ © Mettler-Toledo AG, MarCom, Im Hackacker 15, CH-8902 Urdorf. 03/2009 1
- CH** **Mettler-Toledo (Schweiz) GmbH**, Im Langacher, Postfach, CH-8606 Greifensee
- CN** **Mettler-Toledo Instruments (Shanghai) Co. Ltd.**, 589 Gui Ping Road Cao He Jing,
- D** **Mettler-Toledo GmbH**, Prozeßanalytik, Ockerweg 3, D-35396 Gießen
- F** **Mettler-Toledo, Analyse Industrielle S.A.S.**, 30, Boulevard de Douaumont, BP 949,
- UK** **Mettler-Toledo LTD**, 64 Boston Road Beaumont Leys, GB-Leicester LE4 1AW